

KlentaqLA DNA Polymerase

Amount: 100 μ l (sufficient for 1000 reactions up to 1 kb)

Shipping conditions: Ambient temperature

Storage conditions: -20°C for enzyme, 4°C for 10x Klentaq1 buffer

Thermo stability: Retains at least 85% activity after 1 hour at 95°C

Shelf life: At least 1 year from date of receipt under proper storage conditions.

PRODUCT DESCRIPTION:

KlentaqLA is the Long Accurate version of Klentaq1, a 5'-exonuclease deficient Taq polymerase (an N-terminal deletion of Taq) with improved fidelity and thermostability. 10x buffer composition is: 500 mM Tris-Cl pH 9.2, 160 mM ammonium sulfate, 1% Tween 20, and 35 mM magnesium chloride. We also offer (upon request) 10x buffer at pH 7.9 for better fidelity.

TYPICAL PCR PROTOCOL for a 50 μ l reaction:

Reagent	Volume	Final Concentration
10x Klentaq1 PCR buffer	5 μ l	1x
dNTP mix (10 mM)	1 μ l	200 μ M each
Left Primer	variable	0.2 μ M
Right Primer	variable	0.2 μ M
DNA template [†]	variable	0.1-100 ng
Betaine 5M*	13 μ l (optional)	1.3 M
KlentaqLA Polymerase**	0.1 μ l	2.5 units
de-ionized distilled H ₂ O	Adjust final volume to 50 μ l	-

[†]DNA amount depends mostly on genome size and target gene copy number.

*Betaine is a general PCR enhancer. It usually improves the yield and specificity of amplification especially for longer targets.

**To determine specific optimal enzyme concentration, we strongly recommend an enzyme titration test for each target. Targets larger than 1 kb may require more enzyme.

CYCLING CONDITIONS

1. Denaturing: 94° for 2 minutes for 1 cycle
2. Denaturing: 94° for 40-60 seconds
3. Annealing: 50°-68° depending on the specific T_m primers for 40-60 seconds
4. Extension: 68° for 2 min / 1kb target
5. Repeat steps 2-4 for 25-40 cycles

REFERENCES:

Barnes, W.M. (1994) PCR amplification of up to 35 kb DNA with high fidelity and high yield from bacteriophage templates, PNAS 91, 2216-2220.

U.S. Patent No. 5,436,149

Please visit us on the web at www.klentaq.com for troubleshooting and detailed protocols.

Notice to Purchaser

DNA Polymerase Technology products may not be resold, modified for resale or used to manufacture products without an agreement with DNA Polymerase Technology, Inc. Klentaq1 DNA Polymerase by DNA Polymerase Technology is licensed under US Patent No. 5,436,149. No license for KlentaqLA to be used in a Polymerase Chain Reaction has been purchased by DNA Polymerase Technology, Inc.